



Quick Guide for AutoSwim

1. Download the installation file from the homepage:
www.loligosystems.com/downloads
2. Run the installation program. Follow the instructions on the screen and then restart the PC.
3. Connect the green (wibu) copy protection dongle to a USB port on the PC.
4. Start the AutoSwim program and click *Settings* to verify that all DAQ-BT devices paired with the PC are also listed in AutoSwim.
5. Click *Calibration* to calibrate each swim tunnel/DAQ-BT.
6. Click *Experiment* to start operating the swim tunnel motor.
7. Click *Log* to create a data file and start logging data from a swim tunnel. Please note that each data file logs data from one swim tunnel only.

